




J-KEM® Scientific, Inc.
858 Hodiamont Ave.
St. Louis, MO 63112
(314) 863-5536
Fax (314) 863-6070
E-Mail: jkem911@jkem.com

Precision Pressure Regulator, DPR-1100

J-KEM Scientific's precision vacuum and pressure regulators consist of J-KEM's Infinity-II pressure sensor/controller and KEM-Torr software that presents a graphical user interface to the Infinity controller. The DPR-1100 software displays the system pressure in real-time both graphically and with a digital display, allows the user to construct and run a 12-step pressure ramp, and optionally data logs time, pressure and ramping history.

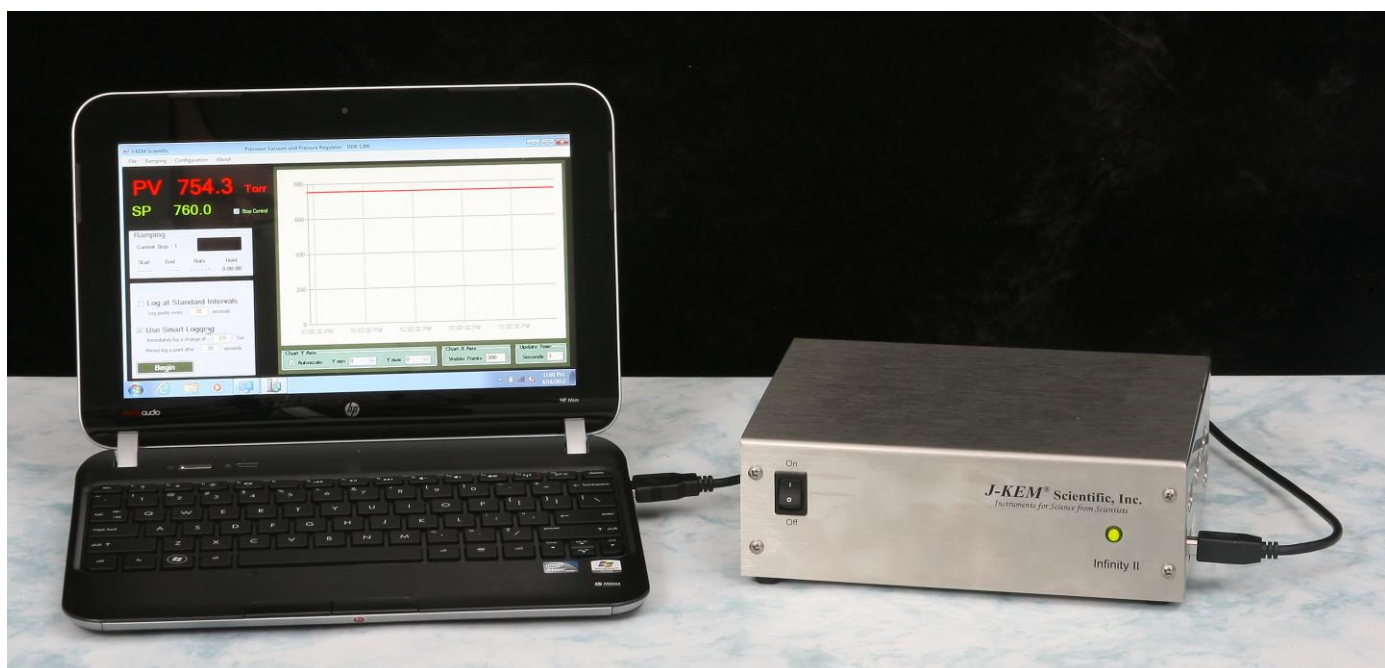


Software Installation

Systems With a PC Controller Pressure regulators ordered with the companion Netbook controller, have KEM-Torr software pre-loaded and no other software setup is required.

System Without a PC Controller KEM-Torr software, PC drivers, and this user manual are included on a flash disk shipped with the Infinity. Do not connect the Infinity controller to your PC until instructed to do so.

1. To install the KEM-Torr software - Insert the flash disk into your PC, open the folder titled Application Software and then double click on the installation icon titled Setup.exe. During the installation, accept all of the default selections.
2. Copy the user manual on the flash disk to the desk top of your PC.
3. Install the Infinity drivers, by opening the folder titled Infinity Drivers and double clicking on the icon titled Infinity_Setup.exe. Allow Windows to install the drivers, and in the process accept all of the default settings.


Hardware Setup for a Pressure Only (single valve) Controller

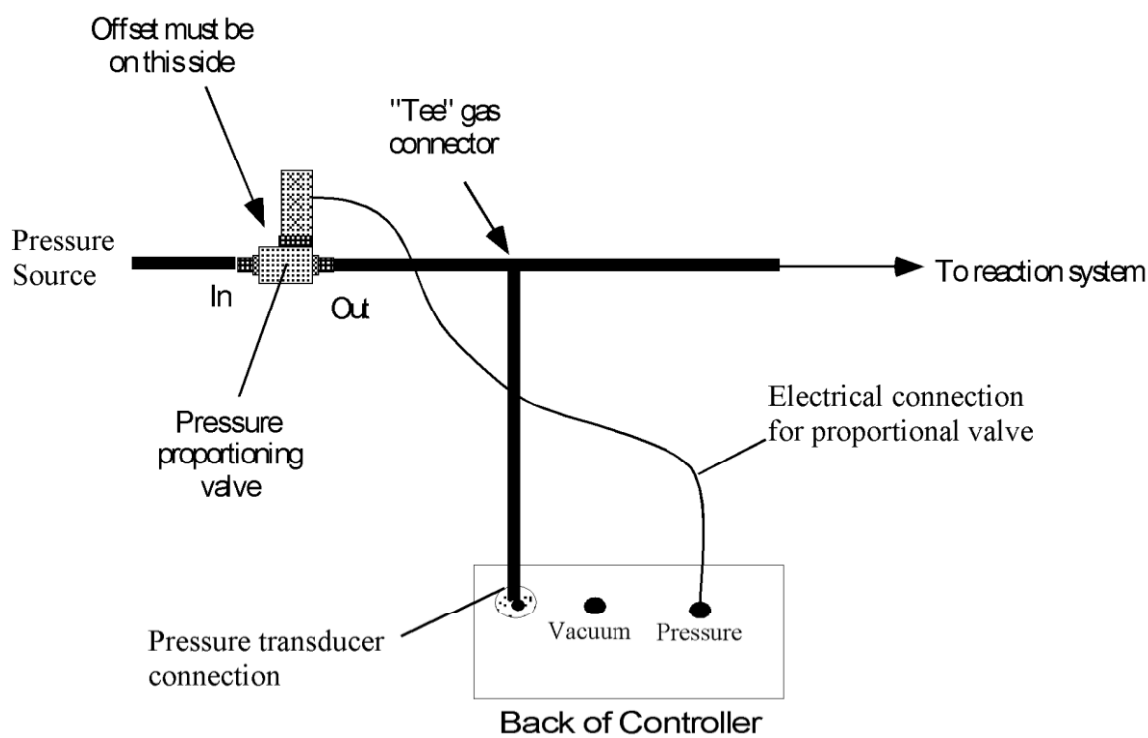

NOTE: The pressure proportioning valve is very sensitive to contaminants that cause the valve to stick. If there is any chance of particulates entering the valve from the reactor, a high efficiency trap must be placed between the proportioning valve and the reaction system to protect the valve. Proportioning valves are not warranted against failure.

1. Connect the inlet of the proportioning valve to the pressure source. The proportioning valve will be one of two types. It will either have a silver coil on top of a silver body (PSV2 - PSV5 valves), or a silver coil on top of a black body (PSV6 – PSV 8 valves). If the valve has a silver body, note that the cylindrical portion of the valve is offset from the square body of the valve, connect the fitting farthest from the cylinder to the pressure source, connect the other fitting to the reaction vessel. If the valve has a black body, connect the port labeled “IN” to the pressure source and the port labeled “Out” to the reactor.
2. Plug the proportioning valve’s gray cord into the electrical connector on the back of the Infinity regulator labeled “Pressure”. Make sure the cable locks into position.
3. Connect the pressure sensing inlet on the back of the Infinity regulator (a threaded or Swagelok fitting) to a Tee connector between the outlet of the valve and the piece of equipment that pressure is being regulated in. Wide bore tubing should be used to make all pressure connections. The length of tubing between the Infinity regulator’s pressure sensing inlet and the reactor under pressure should be kept to a minimum (no longer than 6 feet, but under 3 feet will yield more accurate results). To set up the regulator, simply connect a USB cord between the Infinity-II regulator and the Netbook PC controller or your PC.

Hardware Setup for a Dual Control (dual valve) Controller

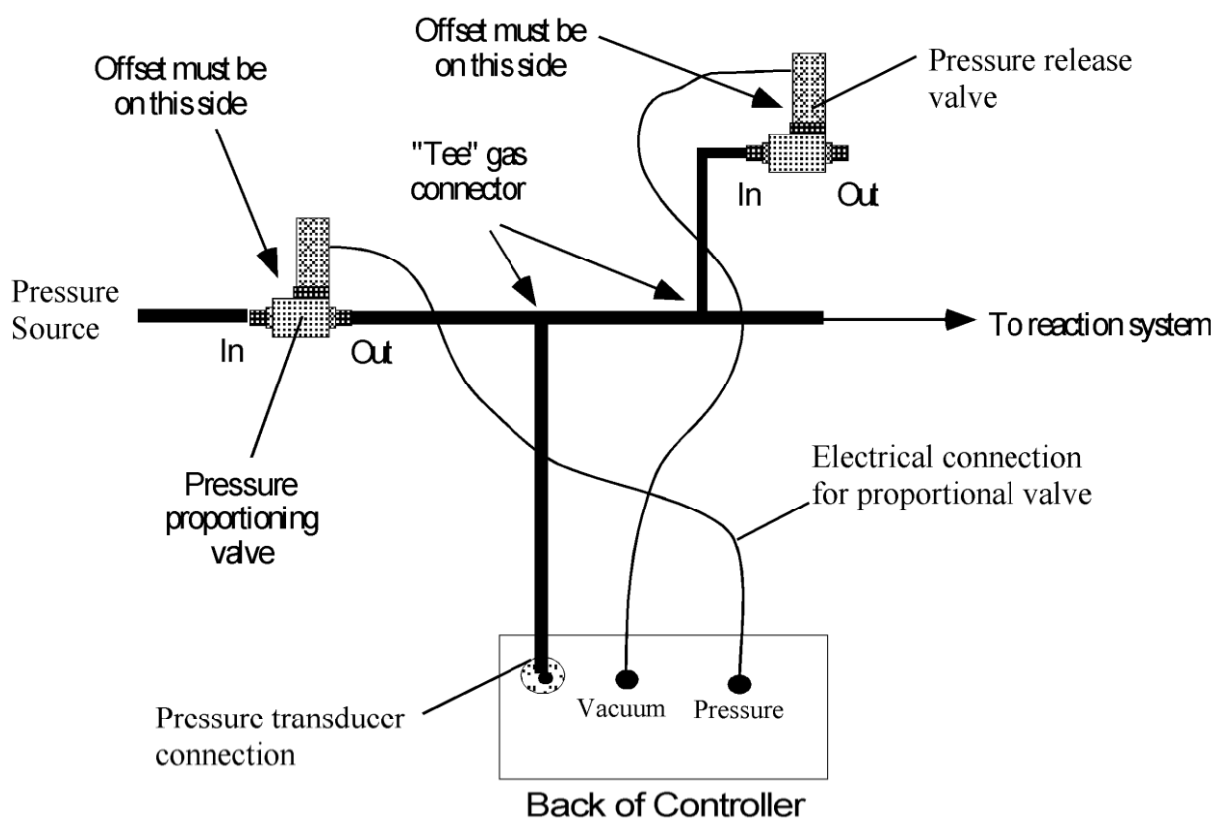


NOTE: The pressurization and pressure release proportioning valves are very sensitive to contaminants which cause the valves to stick. If there is any chance of particulates entering the valve from the reactor, a high efficiency trap must be placed between the proportioning valve and the reaction system to protect the valve. Proportioning valves are not warranted against failure.

1. Connect the inlet of the pressurization proportioning valve to the pressure source. The proportioning valve will be one of two types. It will either have a silver coil on top of a silver body (PSV2 - PSV5 valves), or a silver coil on top of a black body (PSV6 – PSV 8 valves). If the valve has a silver body, note that the cylindrical portion of the valve is offset from the square body of the valve, connect the fitting farthest from the cylinder to the pressure source, connect the other fitting to the reaction vessel. If the valve has a black body, connect the port labeled “IN” to the pressure source and the port labeled “Out” to the reactor.
2. Plug the pressurization proportioning valve’s gray cord into the electrical connector on the back of the Infinity regulator labeled “Pressure”. Make sure the cable locks into position.
3. Connect the inlet of the pressure release proportioning valve to the reactor system using a tee in the inlet pressure line. The proportioning valve will be one of two types. It will either have a silver coil on top of a silver body (PSV2 - PSV5 valves), or a silver coil on top of a black body (PSV6 – PSV 8 valves). If the valve has a silver body, note that the cylindrical portion of the valve is offset from the square body of the valve, connect the fitting farthest from the cylinder to the reactor pressure inlet and leave the other end open to the atmosphere. If the valve has a black body, connect the port labeled “IN” to the reactor pressure inlet and the port labeled “Out” open to the atmosphere.
4. Plug the pressure release proportioning valves gray cord into the electrical connector on the back of the Infinity regulator labeled “Vacuum”. Make sure the cable locks into position.

- Connect the pressure sensing inlet on the back of the Infinity regulator (a threaded or Swagelok fitting) to a Tee connector between the outlet of the valve and the piece of equipment that pressure is being regulated in. Wide bore tubing should be used to make all pressure connections. The length of tubing between the Infinity regulator's pressure sensing inlet and the reactor under pressure should be kept to a minimum (no longer than 6 feet, but under 3 feet will yield more accurate results). To set up the regulator, simply connect a USB cord between the Infinity-II regulator and the Netbook PC controller or your PC.

Software Setup

- Connect a USB cable between the Infinity-II regulator and the Netbook PC controller. Turn on power to the Infinity before starting the KEM-Torr software.
- Start the KEM-Torr software by double clicking on the KEM-Torr icon on the controller's desktop.

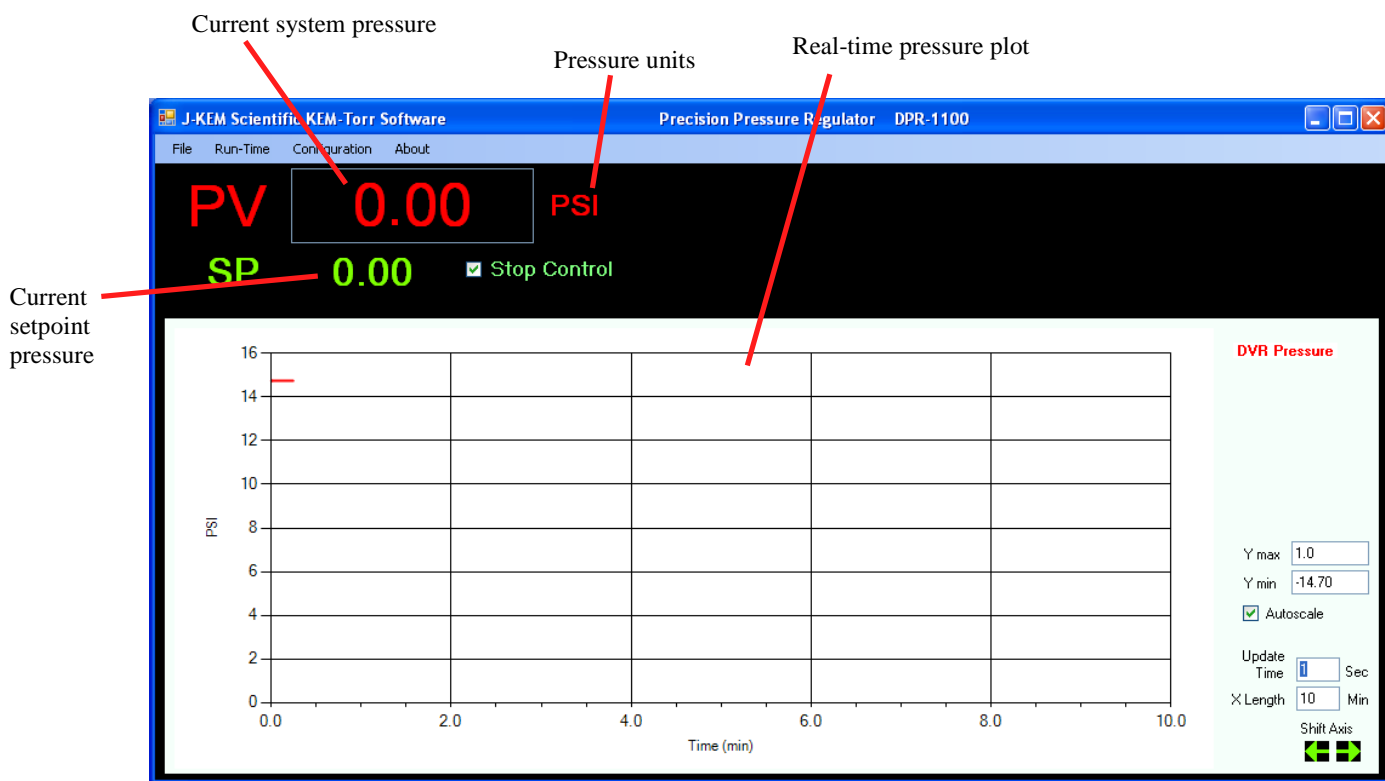

NOTE: The first time the software is started, you may want to set the display configuration of the screen. To set the display units, select Display Units from the Configuration menu. This shows the most common pressure units, but a wider selection of units are available by selecting Customize from the Configuration menu.



The pressure sensed by the regulator is continuously updated and displayed as the PV value in red. Pressures are normally displayed as the absolute pressure in the lab. If the lab's current barometric pressure is below 1.00 atm (760 torr), then a negative pressure is displayed. If the lab's current barometric pressure is above 1.00 atm, then a positive pressure is displayed. If you want the controller to display 0.0 no matter what the lab's current barometric pressure is, click the button titled "Set ATM to 0.0" and the displayed pressure is offset for the lab's current pressure.



Entering a Setpoint Pressure - To enter a setpoint pressure (i.e., the target pressure) place the cursor inside the green setpoint pressure value (i.e., “0.00”) and type the desired setpoint, then hit the Enter (Return) key on the keypad. While the new setpoint is being input, the value turns yellow, when the Return key is hit the new value is entered and the setpoint value turns green again. The regulator begins to operate the pressure proportioning value to adjust the system pressure to the newly entered value.

Single and Dual Valve Controllers – The standard DPR-1100 is programmed to regulate a single pressurization proportioning valve. With only a pressurization valve, the DPR-1100 can increase pressure in the system under test, but it has no ability to decrease pressure. For example, the user can pressurize a chamber to 10 psi and subsequently pressurize it to 12 psi, since the pressure proportioning valve has the ability to increase the system pressure. But if the chamber is pressurized to 10 psi, the DPR-1100 has no ability to decrease the system pressure to 6 psi, since this would require bleeding pressure from the system to the atmosphere. J-KEM does offer an optional feature for the DPR-1100 controller that allows it to both pressurize a system and to depressurize the system. This optional feature requires both a pressurization proportioning valve (the standard valve) and a second depressurization valve (optional). Also the DPR-1100 must be configured for dual valve operation. When the dual valve option is installed, the DPR-1100 can freely increase and decrease pressure in a system under test. Contact J-KEM for information.



Any time during a run, vacuum control can be suspended by placing a check mark in the box titled Stop Control. Removing the check mark restarts vacuum regulation. The system pressure is continuously updated whether the regulator is active or not.

Menu Options – The Run-Time menu contains several options pertaining to an active run.

- Ramping** – Contains the controls needed to create and run a pressure ramp. See the section titled Entering a Vacuum Ramp Program.
- Logging** – Contains controls to configure and initiate logging time and pressure data to a *.csv file on the PC. See the section on Logging for detailed information.
- Charting** – Contains the controls needed to adjust the charting area. See the section titled Charting Controls.
- Fast Update** – When selected, this option provides faster update of the displayed pressure at the cost of slightly more background noise in the pressure signal.

Menu Options – The Configuration menu has several items that affect the operation of the DPR-1100.

- Enable Configuration Menu** – The menu items in the Configuration menu are by default, not enabled to prevent accidentally adjusting a parameter that affect the controller in an adverse manner. To enable the items in the Configuration menu, select Enable Configuration Menu.
- Display Units** – From this menu, the user can select the units that pressure is displayed and entered in. The three common units, PSI, Torr, and Atm are presented in this menu, but other units are available in the Customize menu.
- Monitor DAC Output** – This displays a window on the main form that shows the amount of voltage applied to the proportioning valve. This is for diagnostic purposes only and provides no useable information to the researcher.
- Infinity Configuration** – This opens a form that displays the configuration of the various analog inputs on the Infinity controller. From this form the user can recalibrate the controller if needed, see the section titled Vacuum Regulator Configuration and Calibration.
- Customize** – Allows the user to customize various features of the KEM-Torr software including the units pressures are displayed in, logging defaults and other parameters. See the section titled Customizing the KEM-Torr software.

Entering a Vacuum Ramp Program

Step	Starting Pressure	Ending Pressure	Duration (Min)	Hold Time (Min)	Wait For Setpoint	Step Time
Step 1	760.0	400.0	12.00	0.0	<input type="checkbox"/>	0:12:00
Step 2	500.0	400.0	30.00	0.0	<input checked="" type="checkbox"/>	0:30:00
Step 3	400.0	300.0	30.00	1.0	<input checked="" type="checkbox"/>	0:31:00
Step 4	100.0	50.0	1.00	0.0	<input checked="" type="checkbox"/>	0:01:00
Step 5	50.0				<input type="checkbox"/>	
Step 6					<input type="checkbox"/>	
Step 7					<input type="checkbox"/>	
Step 8					<input type="checkbox"/>	
Step 9					<input type="checkbox"/>	
Step 10					<input type="checkbox"/>	
Step 11					<input type="checkbox"/>	
Step 12					<input type="checkbox"/>	

Total Program Time: **1:14:00**

Save Ramp to Meter

Minimum Pressure: -175.93313486
Maximum Pressure: 175.93313486

Ramp Cycling: Repeat the ramp for cycle(s)

Clear Ramp

The ramping functions of the software support a 12 step vacuum ramp. To display the ramp construction form, select Run-Time -> Ramping -> Build Ramp from the Run-Time menu.

Each row in the table is a single ramp step, a ramp step includes both a pressure change segment and a hold segment.

Starting Pressure – This is the setpoint pressure that the ramp step start at.

Ending Pressure – This is the ending setpoint pressure of the ramp step. For controllers with evacuation control only (single valve controllers), the ending pressure must always be lower than the starting pressure. For evacuation and pressure control units (dual valve controllers), the only restriction of starting and ending pressures is the pressure range of the controller.

Duration (min) – This is the amount of time the program takes to ramp from the Starting to the Ending Pressure.

Hold Time – When a ramp segment completes, the user can optionally enter a hold time for the step. During a hold time, the setpoint is held constant at the ending pressure of the ramp. If you do not want a hold, time, enter a value of 0.

Wait For Setpoint – If you select this option, by checking the checkbox in this column, then the program will not advance to the next ramp step until the sensed pressure in the piece of equipment connected to the DPR-1100 actually reaches the Ending Pressure. For example, if the Starting Pressure is 760 Torr and the Ending Pressure is 400 Torr, the program will perform the instructed ramp step. But, if at the end of the ramp step the pressure in the connected piece of equipment is 420 torr (because the ramp was faster than the piece of equipment could be evacuated), checking this box will cause the ramp program to pause until the pressure in the system actually reaches 400 Torr. When the equipment reaches 400 Torr, then the program continues to the next step. If the box is not checked, then the program continues to the next ramp steps whether the Ending Pressure has been reached or not.

Step Time – This column displays the time required to complete the current ramp step.

When a ramp step completes, including the ramp and hold segments, the ramp program immediately continues at the starting setpoint pressure of the next ramp step. The starting setpoint pressure of the subsequent step is always the ending pressure of the previous step.

When a ramp program completes, the controller remains at the ending setpoint pressure of the last step.

Total Program Time

0:18:30

Save Ramp to Meter

This box shows the total time required to complete all steps of the ramp program. When the ramp program is completely entered, the program is uploaded to the DVR-1000 by clicking the Save Ramp to Meter button.


To delete an entered ramp, you can click the Clear Ramp button on the bottom of the screen.

Ramp Cycling

Repeat the ramp for cycle(s)

If the DVR-1000 is equipped with the optional back-fill pressure option, the Ramp Cycling box is enabled. Ramp Cycling allows the ramp program to restart the ramp program as soon as one pass of the program completes. For example, if a 4-step ramp program is entered, ramp cycling resets the ramp to Step 1 as soon as Step 4 completes. The user can enter then number of times the program will repeat (cycle) in the text box provided.


Storing a Ramp Program - Ramp programs can be stored for later recall. To save a program, select Store Program from the Programs menu. You will be prompted to enter a name for the ramp method. Enter a name, then click OK.

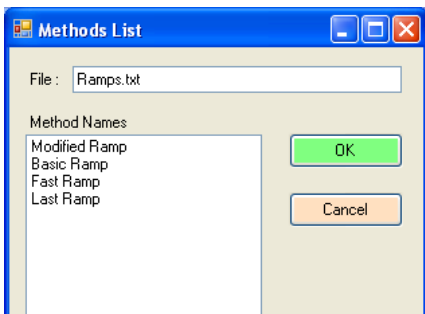

Recalling a Ramp Program – To load a saved ramp, select Recall Program from the Programs menu. A window appears that lists the names of the saved programs, click on the desired program, then click the OK button.

Deleting a Ramp Program – To delete a save program, select Delete a Saved Program from the Programs menu. From the list of programs that appears, click on the program to delete, then click the OK button.

Running a Ramp Program

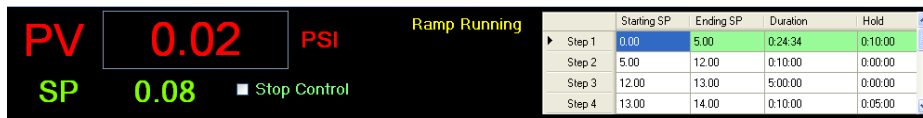

Once a ramp program has been saved to the controller, it appears on the main screen. **Start a Ramp** - To start the ramp, select Ramping -> Start Ramp from the Run-Time menu.

	Starting SP	Ending SP	Duration	Hold
Step 1	0.00	5.00	0:24:08	0:10:00
Step 2	5.00	12.00	0:10:00	0:00:00
Step 3	12.00	13.00	5:00:00	0:00:00
Step 4	13.00	14.00	0:10:00	0:05:00

The current step of a ramp is highlighted in green. During the ramp portion of a program step, the duration time counts down like a digital clock to indicate the time in the ramp remaining. If the program step has a Hold time, then following the completion of the ramp segment, the hold time value begins to count down, indicating the amount of hold time remaining.

	Starting SP	Ending SP	Duration	Hold
Step 1	0.00	5.00	0:23:45	0:10:00
Step 2	5.00	12.00	0:10:00	0:00:00
Step 3	12.00	13.00	5:00:00	0:00:00
Step 4	13.00	14.00	0:10:00	0:05:00


Pause a Ramp - A running ramp can be paused by selecting Ramping -> Pause Ramp from the Run-Time menu. When a ramp is paused, the clock in either the Duration or Hold table cells stop decrementing. To resume the ramp step, select Ramping -> Resume Ramp from the Run-Time menu.

Advance to Next Step – A ramp can be forced to complete the current step and advance to the next step in the program by selecting Ramping -> Advance to Next Step from the Run-Time menu.

Reset Ramp – When a ramp completes, it can be reset and run again by selecting Ramping -> Reset Ramp from the Run-Time menu. When selected, the ramp is reloaded to the meter. To start the ramp, select Ramping -> Start Ramp.

Abort Ramp – A ramp can be terminated by selecting Ramping -> Abort Ramp from the Run-Time menu.

Charting Controls

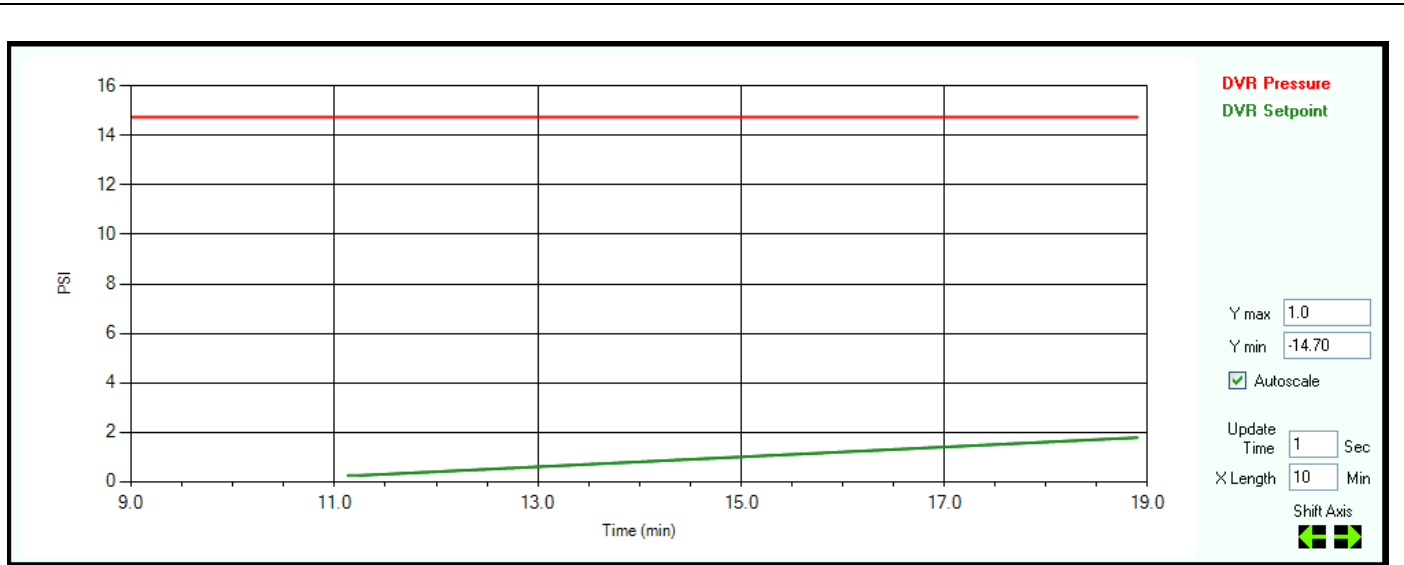

When autoscale is disabled, this sets the minimum value of the Y-axis

When autoscale is disabled, this sets the maximum value of the Y-axis

Allows KEM-Torr to automatically adjust the scale of the Y-axis as needed

Graph update time

Length of the X-axis in units of minutes

Shifts the displayed X-axis left or right

Autoscale - When checked, the graph automatically sets the Y-axis scale to display the plotted data. When unchecked, the user can manually enter the Y-axis minimum and maximum values.

X-Axis Length Sets the length of the X-axis

Update Time Sets the update time of the X-axis and the rate at which data points are collected.


Shift Axis - During normal charting, the axis continuously shifts the X axis so that newly collected points are displayed at the end of the plot. The user can manually shift the displayed plot to view points that have scrolled out of the field of view. Every time the shift left button is clicked, the X plot shifts to older times by the amount specified in the Chart X Axis Length box. Every time the shift right button is clicked, the X plot shifts to newer times by the amount specified in the Chart X Axis Length box. While a plot is shifted left, new points are being logged to the plot, but the plot does not update its view. The view remains static and shows only the section of the chart that has been shifted to. While a charts X-axis is shifted, the back color of the chart changes to light blue.

Menu Options A bit map image of the chart currently displayed can be captured and saved to the PC hard drive. This might be useful for process documentation purposes. To capture the current chart, from the Run-Time menu, select Charting -> Capture Pressure Plot. When this menu option is selected, you are prompted to enter a file name for the plot.

To clear the charting area, select Charting -> Reset. Clearing the charting area only affects the chart, it has not affect on data being logged to a file using the data logging controls.

The setpoint pressure can be plotted on the chart by selecting the option Chart Setpoint Pressure from the Charting menu in the Run-time menu.

Logging Pressure Data

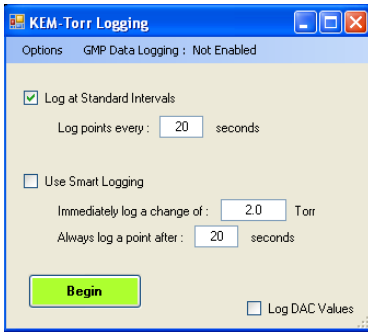




To start logging process data to disk, select Logging, then Start Logging from the Run-Time menu.

There are two logging methods, 1) Standard Logging, and 2) Smart Logging.

Standard Logging is enabled by placing a check mark in the box titled “Log at Standard Intervals”. Standard logging records a data point after every time interval specified in the box titled “Log points every “X” seconds.”

Smart Logging is enabled by placing a check mark in the box titled “Use Smart Logging”. Smart logging records a data point when either of two events occurs 1) the time period set in “Always log a point after “X” seconds” passes, or 2) a change in the process pressure exceeds or equals the value set in “Immediately log a change of “Y” Torr”.


Comparison of Standard and Smart Logging

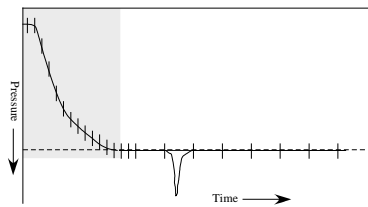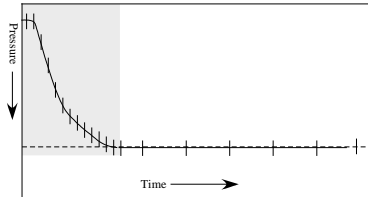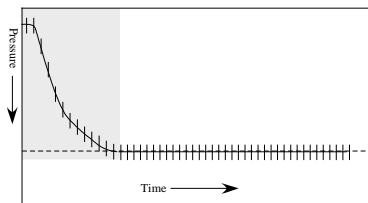

It’s not always necessary, or desirable, to log data points at specific time intervals. For example, consider this plot.


During the initial stage of evacuation it’s desirable to log data points at short, regular intervals (crosses on the plot) in order to define the shape of the curve. But after the pressure stabilizes at the setpoint, rather than logging points at short intervals, which generate large data tables, it’s probably better to log data at longer intervals such as every 1 or 5 minutes as shown in the second plot, because there’s little value in recording a stable pressure reading.

The danger of logging data points at long intervals is that a sharp change in pressure may be missed because it occurred between logged points.

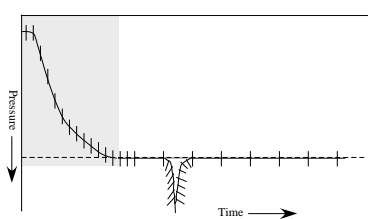

Smart logging addresses this issue by allowing the user to set two parameters. First the user sets the maximum time period between logged points and, second, sets a *log on pressure change value*. If the sensed pressure changes by the *pressure change value* for any reason, this forces a data point to be immediately logged.

With smart logging on, the shape of the pressure spike is perfectly defined because a data point was logged every time the sensed pressure changed by the *pressure change value*. After the sensed pressure stabilizes, data is logged at the longer timer interval. Smart logging provides the best balance between defining the shape of a pressure profile and keeping log files to a manageable size.

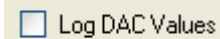


The option to log DAC values is a diagnostic option and is normally not selected. DAC (Digital to Analog Converter) values represent the voltage applied to the vacuum and pressure proportioning valves.

Starting Data Logging

Begin

Click on the Begin button to start data logging. The program prompts the user to enter a file name to save the run data to. The number of logged data points is updated as points are saved.

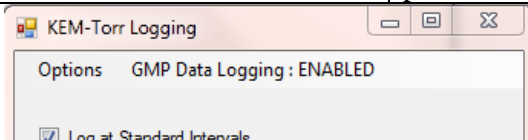


The user has two options for the type of data file that is created. KEM-Torr software can log data to either a compliant or non-compliant data file. Logging data to a GMP compliant data file requires a little more CPU power, but the user will probably not note any change.

Non-GMP Compliant Data File – This is a standard data file that’s logged to Excel in a .csv file format. This file can be open with excel and manipulated in any way.

GMP Compliant Data File – A GMP compliant data file is identical to the non-compliant file, except at the end of the file are appended encrypted hash codes that are unique to each data file. In this case, GMP compliance doesn’t mean that a user is prevented from changing a data point, for example changing a recorded pressure from “760.0 torr” to “300.0 torr”, but that **ANY** change to file is irreversible detectable. KEM-Torr contains a utility that reads the entire data file and then compares that file to the encryption codes saved with the file. If the file matches the encryption codes, then this is a guarantee that no change has occurred to the file since its original creation. If the file does not match the encryption codes, KEM-Net states that the file has been changed since its’ original collection date.

To test a file for data integrity, after the file is closed and no longer logging data, select Run Time -> :Logging -> Log File Integrity -> Verify Log File. You’ll be prompted to select the log file of interest, then KEM-Net will state that the file has not been changed since its original creation, or that it has been changed.


To enable GMP data logging, see the section titled Customizing the DVR-1100 Software.

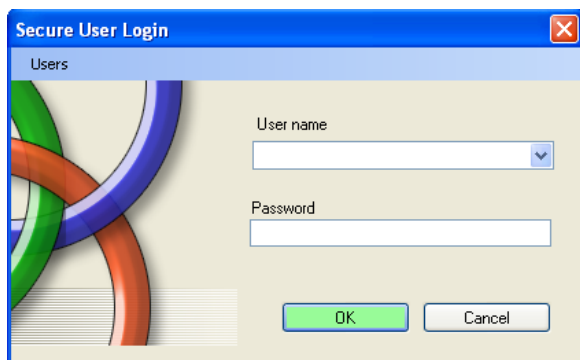

When GMP compliant data logging is selected, the user is prompted to enter their user name and password.


New users can be added by selecting the menu option “Add New Users” from the Users menu. Existing users can be deleted by selecting the menu option “Delete Users” then following the on-screen instructions.

Customizing the DPR-1100 software

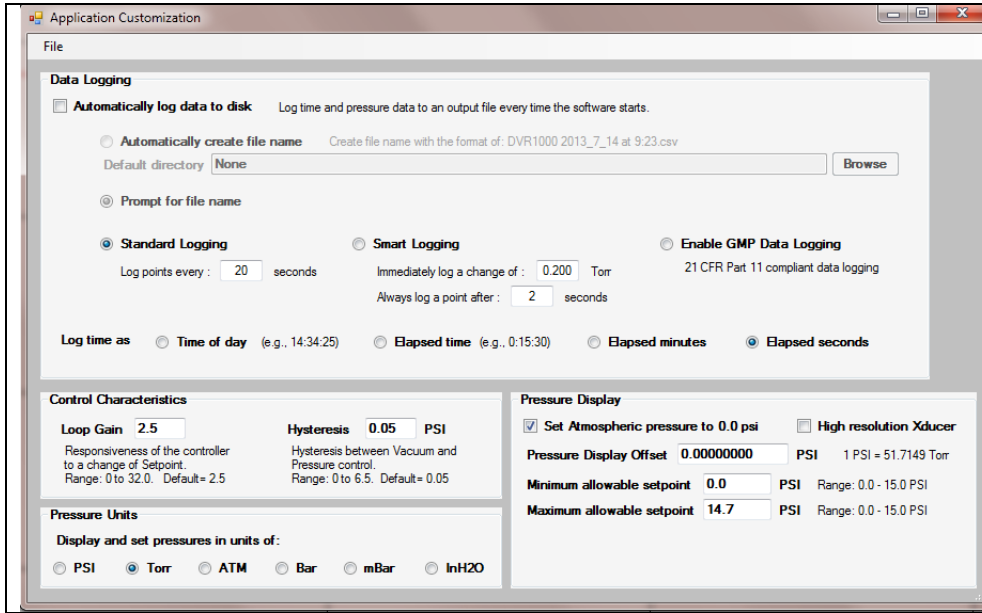


Several parameters of the KEM-Torr can be specified as user defaults. To open the customization screen, select Enable Configuration Menu from the Configuration menu, then select Customize.

Data Logging Controls



<input checked="" type="checkbox"/> Automatically log data to disk	<p>If this option is selected, the controller automatically starts to log time and pressure data to disk as soon as the user starts the control action of the controller by unchecking the “Stop Control”  check box.</p>
<input checked="" type="radio"/> Automatically create file name <input type="radio"/> Prompt for file name	<p>If the option “Automatically create file name” is selected, when data logging starts, the user is not prompted for a file name to save data to, but the system creates a file name based on the current date and time. If this option is not selected, when data logging starts, the user will be prompted for a file name to save data to.</p>
<p>Default Directory</p>	<p>Optionally, the user can enter a default directory to save all data files to.</p>
<p>Log time as:</p>	<p>In the log file, time data can be recorded in any of 4 formats.</p> <p>Time of day – In this format the time the data point is collected is saved at the actual time of day, for example, “4:15 PM”.</p> <p>Elapsed time – This option saves the time that the data point is collected in the format of “Hr : Min : Sec” since the start of the data collection.</p> <p>Elapsed minutes – This option saves the time data point as a decimal value of minutes since the start of the data collection.</p> <p>Elapsed seconds – This option saves the time data point as in integer value of seconds since the start of the data collection.</p>
<p>Type of Data Logging</p>	<p>When logging is set to start automatically, the selection of these variables is used for the logging session.</p> <p>Standard Logging Smart Logging</p> <p>For a discussion on the nature and differences of these two logging methods, see the section above titled Logging Pressure Data.</p>

Enable GMP Data Logging	When data is logged, it can be logged to a simple Excel .csv file in a human-readable, but unsecure file format, or data can be logged into the same file, but with appropriate encryption keys that the file complies with 21 CFR, Part 11 requirements for data logging (GMP Compliant). The file is still human-readable, but appended to the end of the file are codes that will detect if the text of the file is changed in any way. To enable GMP data logging, check the radio button.
--------------------------------	--

Control Characteristics

	The Loop Gain sets how responsive the controller is to a deviation of the
<p>system pressure from the setpoint pressure.. Low values product slow, but stable adjustments to bring the system pressure to the setpoint. High values result in a fast response to a deviation from setpoint, but can also result in unstable control. The default value is 2.5.</p> <p>Hysteresis – Dual valve controllers have a pressure relief and pressurization valve. The algorithm controlling the valves must know how much deviation from the setpoint is required to switch between the two valves. For example, if a flask is being regulated at 6.00 psi, if the flask pressure rose to 6.01 PSI (over pressure), it would not be a good idea to open the pressure relief valve with the expectation that the value would release just the right amount of air to lower the pressure from 6.01 to 6.00 PSI. When the flask is at positive pressures, hysteresis is the amount that the flask pressure must be above the setpoint before the pressure relief valve is activated to correct for the over-pressure condition. The default value for hysteresis is 0.05 PSI. Hysteresis has no affect on how precisely the pressure is regulated in the flask, the Infinity always attempts to regulate the flask pressure to 0.01 PSI.</p>	

Pressure Units

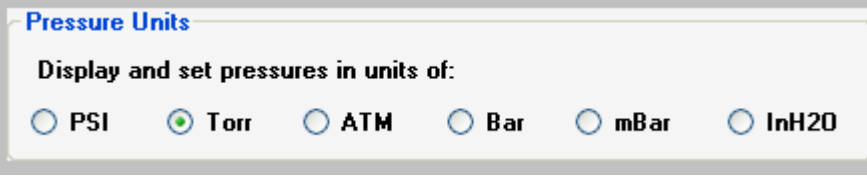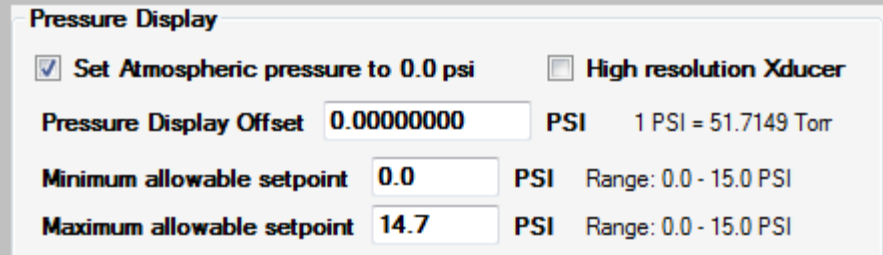
	Specifies the default units for pressure display, data logging, and charting. After selecting the desired unit, the software must be exited and restarted for the change to take effect.
	The user can select to display atmospheric pressure (ATM) as 0.0 or 14.7 psi. If ATM is displayed as 0.0, then system pressures are relative to atmospheric pressure, that is, a setpoint of 1 psi, will be 1 psi above atmospheric pressure, or at 15.7 psi absolute pressure.
<p>If ATM is not displayed as 0.0, then system pressures are displayed as absolute pressures above full vacuum. This option should only be used with DPR-1100's equipped with an absolute style pressure transducer.</p>	

A user offset can be added to the displayed pressure to correct a pressure error present in the display. If a display error is larger than 0.2psi, the controller may need to be recalibrated. Some controllers are supplied with a custom high resolution transducer. Do not check this box unless instructed to do so by J-KEM Scientific.


The user can restrict the pressure setpoint values that can be entered into the controller on the main screen. For example, the user can set the minimum allowable setpoint to a value no lower than 1.5 PSI or to a value no higher than 12.0.

Pressure Regulator Configuration and Calibration


The Infinity board inside of the DPR-1100 is a highly versatile board capable of multiple functions. As can be seen, most of the functionality of the Infinity board is not used for the DPR-1100, but if your application has special requirements, please contact J-KEM for more information.

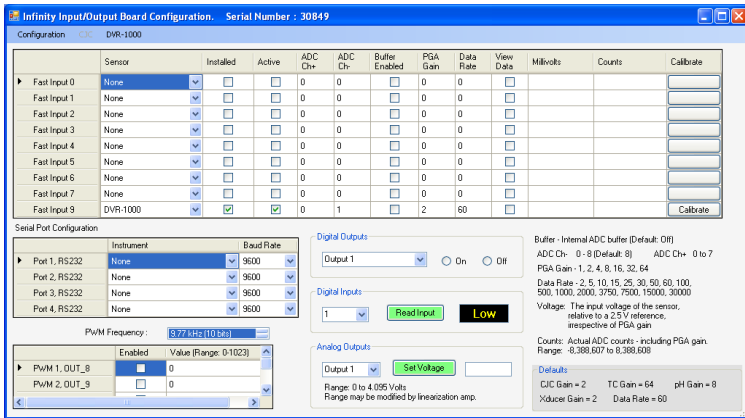

Only one analog input is enabled for the DVR, and that input connects to the pressure transducer. The parameters for this input are locked so that they can not be inadvertently changed. The only control that is not locked is the Calibrate button. Prior to calibrating the controller, you must make sure that you have the appropriate pressure calibrator available.

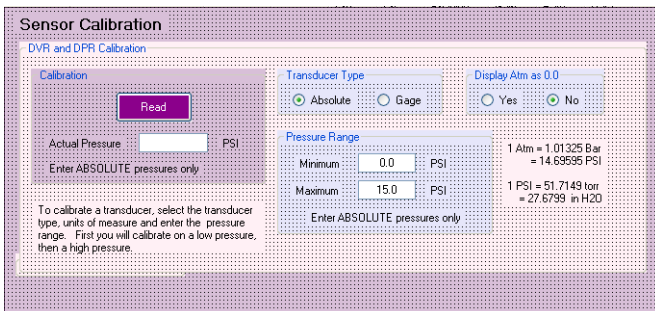

Click the calibration button to open the calibration input screen. In the box titled 'Actual Pressure', enter the value "0.00". After allowing the pressure transducer to sit at atmospheric pressure for 30 seconds, click the Read button to record this pressure point.

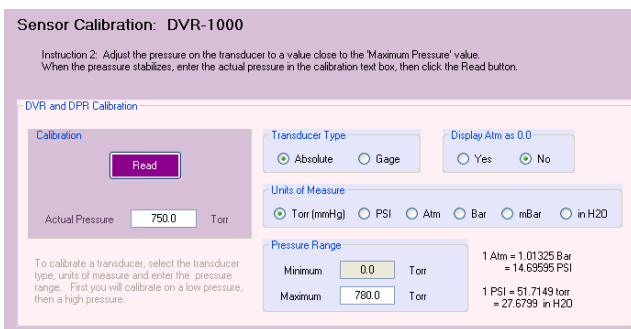

Now connect a stable pressure source to the pressure transducer, but make sure not to exceed the pressure limit of the transducer. Enter the pressure value in the box titled 'Actual Pressure', then click the Read button. The DPR-1100 has a Gage transducer, so enter pressure value as gage pressures.



The regulator now begins to display the sensed pressure on the vacuum transducer. When you are satisfied that the calibrated pressure is reading properly, click the Exit button.

Remote Control of the Infinity Controller

The infinity controller does not require KEM-Torr software to run. For users that want to incorporate the Infinity controller into a larger automation system, the controller can be operated by means of ASCII serial commands.

The controller only accepts one set of communication settings:

Baud: 115200

8 data bits

1 stop bit

No parity

No hardware handshaking.

Understanding the Pressure Unit System Used by the Infinity Controller

Pressure units are entered, and pressure readings are reported in units of “PSI x 10,000”, where atmospheric pressure is reported as 0.00 psi.

So, for example, atmospheric pressure is reported as “0.000”, full vacuum is reported as –146766, and positive 6.5 psi (i.e., 6.5 psi above atmospheric pressure) is reported as 65000.

A pressure of “-53456” is –5.3456 psi, or 483.6 torr $[760.0 - (-5.3456 \text{ psi} * 51.714924 \text{ torr/psi})]$.

Command Set

Note that all commands are case sensitive and are all capital letters. Commands are terminated with carriage return (0x13). The replies to commands, start with the command sent, then any returned data is appended to the returned command string.

Command	Range	Example	Reply	Comments
PCACP	None	PCACP\r	PCACP-1845\r	This command queries the current sensed pressure. The sensed pressure in this example is –1845, or –0.1845 psi. For the DVR-1000 and DDR-1200, pressures are reported as Absolute pressures, while the DPR-1100 reports pressures as gage pressures.
PCASP	-150,000 to 2,500,00	PCASP-100000\r	PCASP-100000\r	This command both sets or queries the current setpoint pressure. Pressure units are ABSOLUTE pressure in PSI x 10,000. When no value follows the PCASP command, the current setpoint pressure is returned in the reply. When a value follows the PCASP command, this sets the current setpoint pressure to the passed value. A command of PCASP-100000 sets the setpoint of –10.00 psi , or 242.9 torr.
	None	PCASP\r	PCASP-10034\r	Queries the current setpoint pressure.

PCARN	0 to 3	PCARN2\r	PCARN2\r	This command starts or stops pressure control. The argument passed with the command depends on the type of Infinity controller in use. Digital Vacuum Regulator, DVR-1000 – If the regulator is fitted only with a vacuum proportioning valve, the command PCARN2 starts vacuum control. If the DVR-1000 has the optional back-fill valve, then the command PCARN3 starts vacuum control and back-fill control. Digital Pressure Regulator, DPR-1100 – The command PCARN1 starts pressure regulation. Dual Range Pressure Regulator, DDR-1200 – The command PCARN3 starts both vacuum and pressure regulation. The command PCARN0 stops pressure regulation in all model numbers.
PWM4D	None	PWM4D\r	PWM4D692\r	This returns a digital value that is representative of the voltage applied to the vacuum proportioning valve. A value of 0 means no voltage is applied to the valve, a value of 4095 means that 24 Vdc is applied to the valve.
PWM2D	None	PWM2D\r	PWM2D692\r	This returns a digital value that is representative of the voltage applied to the pressure proportioning valve. A value of 0 means no voltage is applied to the valve, a value of 4095 means that 24 Vdc is applied to the valve.
PINFO	None	PINFO\r	MFC V3.1,12-Apr-13,4.116\r	Queries product information and program date. This is a good command to verify that you have serial communication with the controller
PCAHP	-150,000 to 2,500,000	PCAHP-140987\r	PCAHP-140987\r	Set the low pressure calibration point. To calibrate the controller, 1) introduce a known low pressure point, (close to full vacuum) and then send the pressure in PSI x 10,000 as an absolute pressure. So, for example, if the low pressure standard point is 1.2 torr, then send the command: PCAHP-146727\r
PCALP	-150,000 to 2,500,000	PCALP-213\r	PCALP-213\r	Set the high pressure calibration point. To calibrate the controller, 1) introduce a known pressure point close to atmospheric pressure and then send the pressure in PSI x 10,000 as an absolute pressure. So, for example, if the high pressure standard point is 758.9 torr, then send the command: PCAHP-213\r

Here is a reasonable startup sequence:

PINFO - Tests to make sure you have serial communications

PCASP##### - Send in a new setpoint

PCARN3 (or PCARN1, or PCARN2. Use the appropriate command) - Start pressure control

PCACP - The command queries the current sensed pressure.

New pressure setpoints can be sent at any time

To create a ramp, send in a timed sequence of setpoints to create the desired pressure ramp.

To stop pressure control and close both the vacuum and pressure control valves, send the command PCARN0.